

*n*Taq-Pure (purified from insect cell)

Cat.#	Size	Conc.
P025PI	250 units	5 units/μl
P050PI	500 units	5 units/μl

Store at -20°C

Supplied with: 10X *n*Taq Buffer (Mg
dNTP Mixture (2 mM each)
Sterile water

India Contact:

Life Technologies (India) Pvt. Ltd.
306, Aggarwal City Mall, Opposite M2K Pitampura,
Delhi - 110034 (INDIA).
Ph: +91-11-42208000, 42208111, 42208222
Mobile: +91-9810521400
Fax: +91-11-42208444
Email: customerservice@lifetechindia.com
Web: www.lifetechindia.com

Product description

Clean *n*Taq DNA polymerase is a highly purified polymerase especially for PCR reactions where freedom from bacterial genomic DNA is essential. A gene encoding the *Thermus aquaticus* (Taq) DNA polymerase was cloned and expressed in *Spodoptera frugiperda* (Armyworm), and the enzyme was purified to homogeneity. The purified Taq polymerase (*n*Taq) has optimal activity at high temperatures (72°C), which helps amplify secondary-structured regions. *n*Taq DNA Polymerase has greatly reduced activity of 5'→3' exonuclease activity. As a result, formation of erroneously amplified products caused by the 5'→3' exonuclease activity of wild type polymerase is effectively prevented. In addition, because this enzyme was purified from completely different organism to bacteria (such as *E. coli*), no prokaryotic DNA contamination exist.

Characteristics

- Absolutely free from *E. coli* genomic DNA contamination
- Enhanced sensitivity
- Molecular weight: 94 kDa
- Error rate: 2.4×10^{-5}
- Thermal stability: Half life of 40 min at 95°C
- A-tail formation at 3' ends of amplified duplex DNA

10X *n*Taq Buffer (Mg²⁺ plus)

containing 15 mM MgCl₂

For Research Use Only. Not for use in diagnostic procedures.

ISO9001 ISO14001 ISO13485

*n*Taq-Pure (purified from insect cell)

Cat.#	Size	Conc.
P025PI	250 units	5 units/μl
P050PI	500 units	5 units/μl

Store at -20°C

Supplied with: 10X *n*Taq Buffer (Mg²⁺ plus)
dNTP Mixture (2 mM each)
Sterile water

India Contact:

Life Technologies (India) Pvt. Ltd.
306, Aggarwal City Mall, Opposite M2K Pitampura,
Delhi - 110034 (INDIA).
Ph: +91-11-42208000, 42208111, 42208222
Mobile: +91-9810521400
Fax: +91-11-42208444
Email: customerservice@lifetechindia.com
Web: www.lifetechindia.com

Product description

Clean *n*Taq DNA polymerase is a highly purified polymerase especially for PCR reactions where freedom from bacterial genomic DNA is essential. A gene encoding the *Thermus aquaticus* (Taq) DNA polymerase was cloned and expressed in *Spodoptera frugiperda* (Armyworm), and the enzyme was purified to homogeneity. The purified Taq polymerase (*n*Taq) has optimal activity at high temperatures (72°C), which helps amplify secondary-structured regions. *n*Taq DNA Polymerase has greatly reduced activity of 5'→3' exonuclease activity. As a result, formation of erroneously amplified products caused by the 5'→3' exonuclease activity of wild type polymerase is effectively prevented. In addition, because this enzyme was purified from completely different organism to bacteria (such as *E. coli*), no prokaryotic DNA contamination exist.

Characteristics

- Absolutely free from *E. coli* genomic DNA contamination
- Enhanced sensitivity
- Molecular weight: 94 kDa
- Error rate: 2.4×10^{-5}
- Thermal stability: Half life of 40 min at 95°C
- A-tail formation at 3' ends of amplified duplex DNA

10X *n*Taq Buffer (Mg²⁺ plus)

containing 15 mM MgCl₂

For Research Use Only. Not for use in diagnostic procedures.

ISO9001 ISO14001 ISO13485

Your Molecular & Cell Technology Partner

Standard PCR conditions

- PCR mixture^a

10X <i>n</i> Taq Buffer (Mg ²⁺ plus)	2 μl
<i>n</i> Taq-Pure DNA Polymerase ^b (5 units/μl)	0.2 μl
dNTP Mixture (2 mM each)	2 μl
Template DNA ^c (0.1~500 ng/μl)	1 μl
Primer 1 (5 pmole/μl)	1 μl
Primer 2 (5 pmole/μl)	1 μl
Sterile water	up to 20 μl
^a Assemble the reaction mixture on ice	
^b Add the PCR polymerase at the final step	
^c Plasmid DNA, 0.1 ng~30 ng; genomic DNA, 50 ng~500 ng	

- PCR cycle

Initial denaturation ^a	95°C	2 min
Denaturation	95°C	30 sec
Annealing ^b	55°C~65°C	30~60 sec
Elongation	72°C	1 min/kb
Number of cycles	25~35 times	
Final elongation	72°C	5 min
^a When cycles are over, keep the reaction mixture at 4°C; may add 10 mM EDTA until use to prevent DNA degradation.		
^b Recommended annealing temperatures is 5 to 10°C below the lower T _m of the two primers used.		

Applications

- Standard / General PCR
- PCR with bacterial DNA
- *E. coli* contamination studies
- Microbial (i.e. 16S/23S) screening studies
- Forensic studies
- PCR cloning
- RT-PCR

Quality control

- Purity: >99% on SDS-PAGE
- Endonuclease-free
- Exonuclease-free
- RNase-free
- Inhibitor-free

Unit definition

One unit is defined as the amount of enzyme required to incorporate 10 nmol of dNTP into acid insoluble materials in 30 min at 74°C in a 50 μl reaction mixture (20 mM Tris-HCl/ pH 8.8, 50 mM KCl, 2.5 mM MgCl₂, 10 mM β-mercaptoethanol, 12.5 μg of calf thymus DNA).

Storage buffer

20 mM Tris-HCl (pH 7.9), 100 mM KCl, 0.1 mM EDTA, 1 mM DTT, 0.5% NP-40, 0.5% Tween-20, 50% glycerol.

Applications

- Standard / General PCR
- PCR with bacterial DNA
- *E. coli* contamination studies
- Microbial (i.e. 16S/23S) screening studies
- Forensic studies
- PCR cloning
- RT-PCR

Quality control

- Purity: >99% on SDS-PAGE
- Endonuclease-free
- Exonuclease-free
- RNase-free
- Inhibitor-free

Unit definition

One unit is defined as the amount of enzyme required to incorporate 10 nmol of dNTP into acid insoluble materials in 30 min at 74°C in a 50 μl reaction mixture (20 mM Tris-HCl/ pH 8.8, 50 mM KCl, 2.5 mM MgCl₂, 10 mM β-mercaptoethanol, 12.5 μg of calf thymus DNA).

Storage buffer

20 mM Tris-HCl (pH 7.9), 100 mM KCl, 0.1 mM EDTA, 1 mM DTT, 0.5% NP-40, 0.5% Tween-20, 50% glycerol.

Your Molecular & Cell Technology Partner

Standard PCR conditions

- PCR mixture^a

10X <i>n</i> Taq Buffer (Mg ²⁺ plus)	2 μl
<i>n</i> Taq-Pure DNA Polymerase ^b (5 units/μl)	0.2 μl
dNTP Mixture (2 mM each)	2 μl
Template DNA ^c (0.1~500 ng/μl)	1 μl
Primer 1 (5 pmole/μl)	1 μl
Primer 2 (5 pmole/μl)	1 μl
Sterile water	up to 20 μl
^a Assemble the reaction mixture on ice	
^b Add the PCR polymerase at the final step	
^c Plasmid DNA, 0.1 ng~30 ng; genomic DNA, 50 ng~500 ng	

- PCR cycle

Initial denaturation ^a	95°C	2 min
Denaturation	95°C	30 sec
Annealing ^b	55°C~65°C	30~60 sec
Elongation	72°C	1 min/kb
Number of cycles	25~35 times	
Final elongation	72°C	5 min
^a When cycles are over, keep the reaction mixture at 4°C; may add 10 mM EDTA until use to prevent DNA degradation.		
^b Recommended annealing temperatures is 5 to 10°C below the lower T _m of the two primers used.		